



# Product datasheet FERMACIDAL D2

Disinfectant for surfaces, objects, tools - Solution ready to use

**Ingredients:** Quarternary ammonium compounds. Does not contain alcohol, mercury, formaldehyde, phenol or other toxic compounds.

**Active agents:**

Alcyl(C12-18)dimethylbenzylammoniumchloride  
(ADBAC) – 0,67 mg/ml

Alcyl(C12-14)dimethyl(ethylbenzyl)ammoniumchloride  
(ADEBAC) – 0,67 mg/ml

Didecyldimethylammoniumchloride  
(DDAC) – 0,67 mg/ml

**Properties:** cleaning, blood dissolving, neutralising bacterial odour, does not spot, suitable for use on most surface materials, non-corrosive, odourless, non-irritant, non-volatile, skin-friendly, non-toxic for humans and animals as well as for mammalian cells and tissues in culture, biodegradable.

**Efficiency:** Bactericidal (E. coli), fungicidal (Candida), efficient against mycobacteria (TbB), mycoplasma, hepatitis B, HIV, and other viruses (e.g. rota, influenza A, H5N1, H1N1, corona, etc.), prevents endospores from flourishing.

**Safety information:** Disposal of the biocidal product and its packaging according to official regulations.  
P102 – Keep out of reach of children. Harmful if not used as directed.

**Fields of application:** For contamination prevention in cell culture- and molecularbiology laboratories, hospitals, therapy- and nursing facilities, retirement homes, schools, kindergartens, solariums, pedicure- and cosmetic studios, car rentals, restaurants, hotels, canteens, shops, food processing industry, public buildings (bathhouses, sports facilities, camping, etc.).

Surfaces, objects and tools (cell and tissue culture incubators, clean benches, furniture, devices, instruments, tools, tables, handles, handrails, tapware, etc.)

**Instructions for use:** spray objects daily until wet. Then let surfaces air-dry. Drying time = Reaction time, no wiping!

**Registered as biocide:**

**BAG / FOPH No. CHZB0445**

**BauA Reg.-No. N 24040**

**VAH-listed**

**Manufacturer:** IC Products SA, CH-6596 Gordola

